



Benefit of Using Yamazen Inject Column Drastic Column Downsizing Is Possible.

- If a small separation column is used with an Inject column, this can purify a sample as well as a large separation column used alone.
- A drastic reduction in solvent use


Sample: Butyl-P-hydroxybenzoate (200mg)
Methyl-P-hydroxybenzoate (200mg)

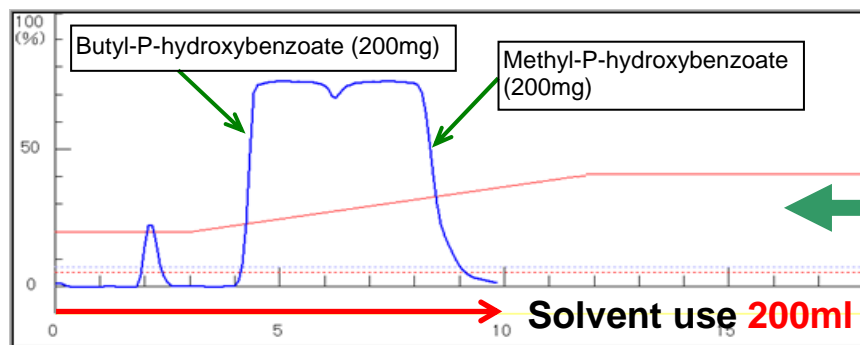

Only an Universal column L (40g) is used.

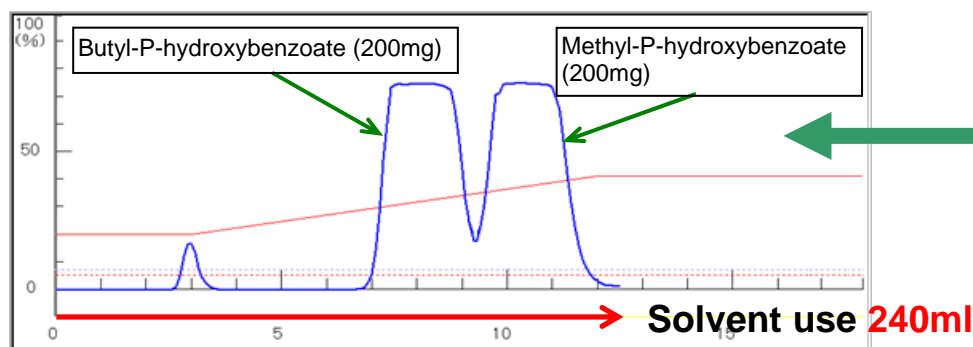

A pair of an Universal column L (40g) and an Inject column M for sample loading are used.

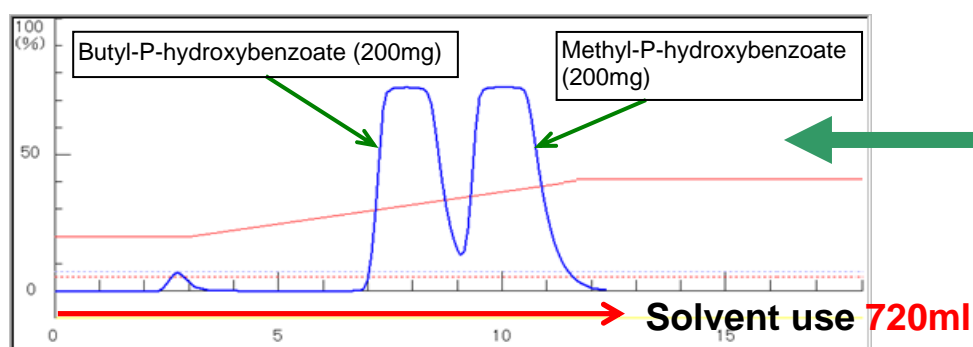

Only an Universal column 3L (133g) is used.

A pair of an Universal column L (40g) and a M size Inject column can purify a sample as well as an Universal column 3L (133g) is used alone.